

Energy Transfer in a Liquid Scintillator Investigated using Time-Resolved X-ray Excited Luminescence Spectroscopy

AN_P71; Stuart Thomson

Introduction

Liquid scintillator detectors are widely utilised in neutrino and astroparticle detection experiments. One liquid scintillator that is increasingly being used in large scale detection experiments is linear alkylbenzene (LAB). LAB is an attractive scintillator due to its low cost, high flash point and low toxicity; making it easier to handle than previously used toxic and flammable organic solvents. LAB is commonly used together with the fluor 2,5-diphenyloxazole (PPO) which increases luminescence yield and shifts the luminescence to longer wavelengths. The LAB/PPO scintillator system is already used at the Daya-Bay¹ and RENO² neutrino detection experiments and is the scintillator of choice for the upcoming SNO+³ and JUNO⁴ detectors.

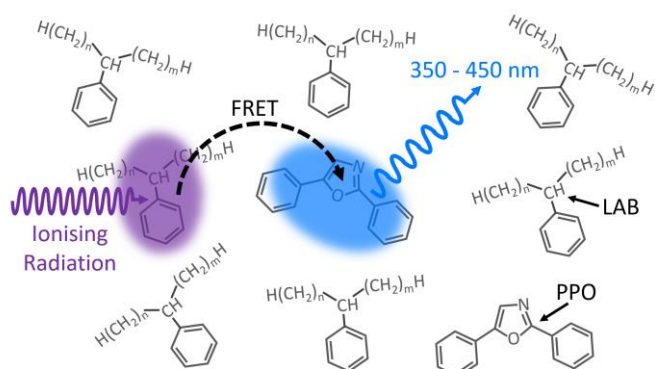


Figure 1: LAB/PPO Scintillator System.

A schematic of the scintillation process in the LAB/PPO system is shown in Figure 1. Ionising radiation, such as γ -rays, travel through the LAB solvent and excite a molecule of LAB to a high energy state. The excess energy in the LAB is non-radiatively transferred (FRET) to a nearby PPO molecule promoting it to an excited state. Radiative de-excitation of the PPO back to its ground state results in violet/blue luminescence which can be detected.

Since energy is transferred between the LAB and PPO via FRET, the luminescence behaviour will depend on the number of PPO molecules and their distance from excited LAB molecules. Higher concentrations of PPO decrease the average spacing between excited LAB and PPO molecules, increasing the luminescence yield, and decreasing the average luminescence lifetime. Many detection experiments utilise pulse shape discrimination techniques, which analyse the luminescence decay profile after a scintillation event, to distinguish different radioactive particles.

Understanding the luminescence decay behaviour of the scintillator system is crucial for this type of analysis. In this application note, the effect of PPO concentration on the luminescence decay profile of the LAB/PPO system is investigated using time-resolved X-ray excited luminescence spectroscopy with an Edinburgh Instruments FLS1000 Photoluminescence Spectrometer.

Materials and Methods

PPO (Sigma Aldrich) was dissolved in LAB (Cepsa Canada) at a series of concentrations between 1 g/L and 20 g/L, and pipetted into quartz cuvettes. X-ray excited luminescence spectra and decays of the LAB/PPO solutions were measured using an Edinburgh Instruments FLS1000 Photoluminescence Spectrometer coupled to the XS1 X-ray Sample Chamber (Figure 2).

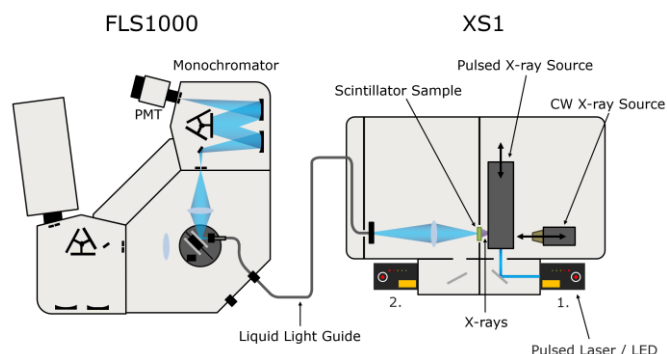


Figure 2: FLS1000 & XS1 X-ray excited luminescence spectroscopy setup.

The XS1 X-ray Sample Chamber is an X-ray shielded enclosure that can be equipped with multiple X-ray sources for the investigation of new scintillator materials. To acquire the X-ray excited luminescence spectra of LAB/PPO the XS1 was equipped with a compact 60 kV continuous-wave X-ray source. For the luminescence decay measurements, the LAB/PPO was excited by a 40 kV pulsed X-ray source which was optically pumped using an Edinburgh Instruments HPL-450 pulsed diode laser. The X-ray excited luminescence from the scintillator is collected by a liquid light guide which delivers it to the FLS1000 where it is wavelength selected and detected. The FLS1000 was equipped with a PMT-900 detector that is sensitive from 200-870 nm and time-correlated single-photon counting (TCSPC) electronics.

Energy Transfer in a Liquid Scintillator Investigated using Time-Resolved X-ray Excited Luminescence Spectroscopy

AN_P71; Stuart Thomson

Results and discussion

The X-ray excited luminescence spectrum of LAB/PPO was first acquired using the CW X-ray source and is shown in Figure 3. The luminescence has a peak at 365 nm and originates purely from the PPO component with no luminescence directly from LAB.

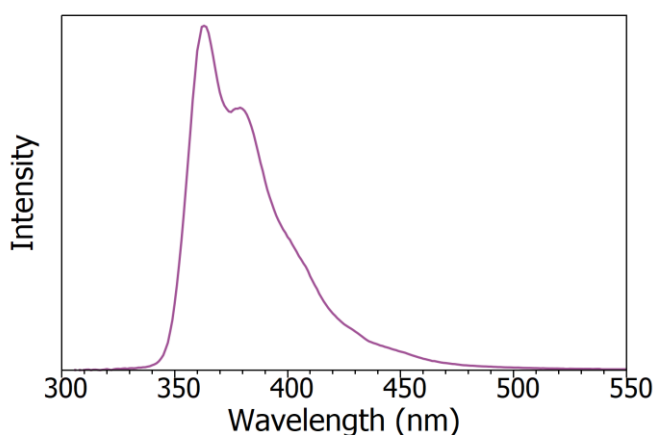


Figure 3: X-ray excited luminescence spectrum of LAB/PPO (8 g/L).

Next, the effect of PPO concentration on the luminescence decay profile of LAB/PPO was investigated. Eight LAB/PPO solutions with varying concentrations of PPO were excited using the 40 kV pulsed X-ray source and their decay measured using TCSPC (Figure 4). The shape of the decay profiles is complex, comprising of an initial prompt component followed by a long tail. The prompt component can be assigned to radiative de-excitation of PPO from its singlet excited state to the singlet ground state which is an allowed transition and therefore occurs rapidly. In contrast, molecules that have been excited into their triplet states cannot de-excite directly to the ground state and instead proceed through alternative de-excitation channels, such as triplet-triplet annihilation, which gives rise to the long tail in the luminescence decays.

To extract the decay components the luminescence decays were fit with either a two (1 - 6 g/L) or three (8 - 20 g/L) exponential decay model using reconvolution fitting with Edinburgh Instruments FAST lifetime analysis software (solid lines in Figure 4). The lifetime and intensity fraction of each decay component is given in Table 1. The fastest decay component (τ_1) corresponds to the radiative singlet de-excitation of PPO and is the dominant contribution to the total luminescence for all PPO concentrations, with a luminescence fraction between 72 and 96 %.

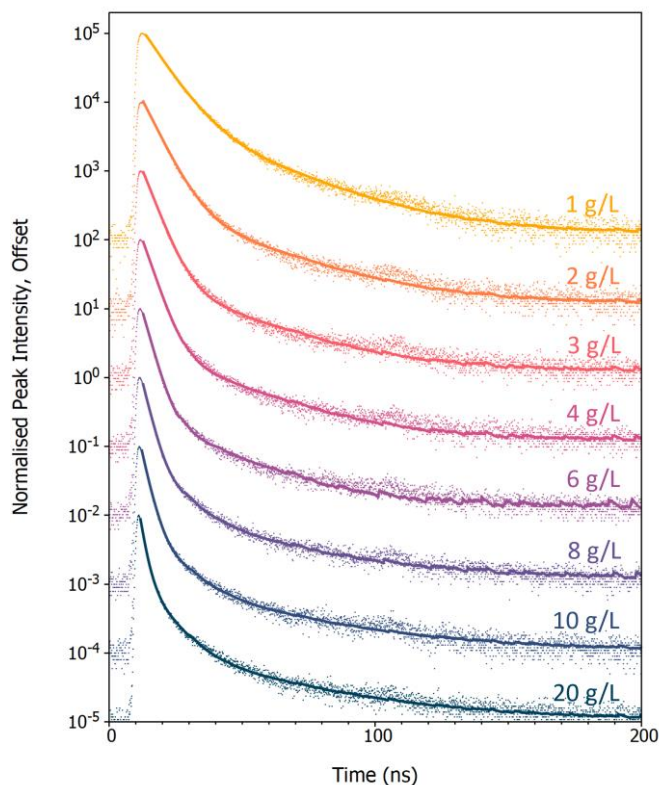


Figure 4: X-ray excited luminescence decays of LAB/PPO at varying PPO concentrations acquired using TCSPC. The scatter plots are the experimental TCSPC decays and solid lines are a two or three exponential component best fit to each decay.

Table 1: Decay times (τ_i) and decay component fractions (f_i) for varying PPO concentrations in LAB after pulsed X-ray excitation.

PPO Conc.	τ_1 (ns)	f_1 (%)	τ_2 (ns)	f_2 (%)	τ_3 (ns)	f_3 (%)
1 g/L	7.1	85.8	26.6	14.2		
2 g/L	5.4	90.9	28.3	9.1		
3 g/L	4.3	94.1	27.2	4.3		
4 g/L	3.7	94.2	26.5	5.8		
6 g/L	3.0	96.2	22.9	3.8		
8 g/L	2.5	89.5	9.3	7.1	36.9	3.4
10 g/L	2.3	79.9	8.8	13.0	39.7	7.1
20 g/L	1.7	71.8	7.5	19.9	38.5	8.3

Energy Transfer in a Liquid Scintillator Investigated using Time-Resolved X-ray Excited Luminescence Spectroscopy



AN_P71; Stuart Thomson

Figure 5 shows the evolution of τ_1 with PPO concentration. Since the excitation of a PPO molecule proceeds via non-radiative energy transfer from an excited LAB molecule τ_1 is dependent on the energy transfer rate. As the concentration of PPO is increased the average spacing between an excited LAB molecule and a neighbouring PPO molecule decreases which increases the energy transfer rate resulting in the decrease in τ_1 with concentration seen in Figure 5. At the highest PPO concentrations, energy transfer is no longer the rate limiting step and τ_1 approaches the intrinsic radiative lifetime of PPO in LAB which is < 2 ns.⁶

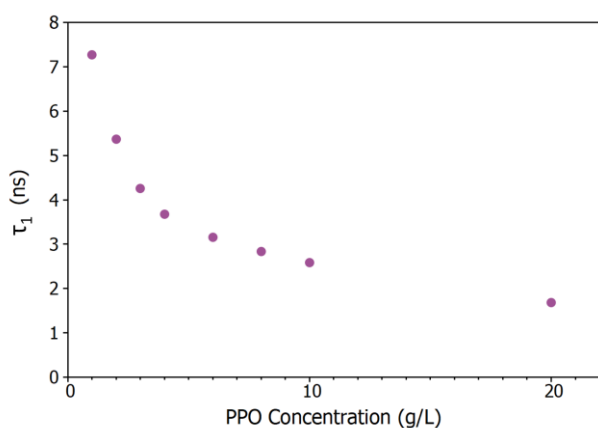


Figure 5: Variation of τ_1 in LAB/PPO with PPO concentration.

Conclusion

The X-ray excited luminescence decay properties of LAB/PPO at varying PPO concentrations were investigated using an Edinburgh Instruments FLS1000 with the XS1 X-ray Sample Chamber Accessory. The lifetime of the primary decay component was found to decrease from 7.1 ns to 1.7 ns as the

concentration of PPO was increased from 1 to 20 g/L. The XS1 X-ray Sample Chamber accessory extends the functionality of the FLS1000 PL spectrometer into the X-ray region, creating a powerful characterisation tool for the development of new scintillator materials.

Acknowledgements

We are grateful to Dr. Sasmit Gokhale from Brookhaven National Laboratory for providing the LAB and PPO used in this application note.

References

1. F. An et al., The Daya Bay Collaboration, Observation of Electron-Antineutrino Disappearance at Daya Bay, *Phys. Rev. Lett.* **108** 171803 (2012)
2. J. Ahn et al., The RENO Collaboration, Observation of Reactor Electron Antineutrinos Disappearance in the RENO Experiment, *Phys. Rev. Lett.* **108** 191802 (2012)
3. S. Andringa et al., The SNO+ Collaboration, Current Status and Future Prospects of the SNO+ Experiment, *Adv. High Energy Phys.* **2016** 6194250 (2016)
4. F. An et al., The JUNO Collaboration, Neutrino Physics with JUNO, *J. Phys. G* **43** 030401 (2016)
5. H. M. O'Keffe et al. Scintillation Decay Time and Pulse Shape Discrimination in Oxygenated and Deoxygenated Solutions of Linear Alkylbenzene for the SNO+ Experiment, *Nucl. Instrum. Methods Phys. Res. A* **640** 119-122 (2011)
6. T. M. Undagoitia et al., Fluorescence Decay-time Constants in Organic Liquid Scintillators, *Rev. Sci. Instr.* **80** 043301 (2009)



For more information, please contact:

+44 (0) 1506 425 300
sales@edinst.com
www.edinst.com