### High-Resolution Raman & PL Imaging of MoS<sub>2</sub>

AN\_R15; Angela Flack



#### Introduction

Transition metal dichalcogenides (TMDs) are a class of 2D materials which have unique optical and electronic properties. They are semiconductors with the chemical structure  $MX_2$  where M is a transition metal and X a chalcogen, Figure 1. After the isolation of graphene layers, research into alternative 2D layered materials which offer the same, or better, properties has surged.

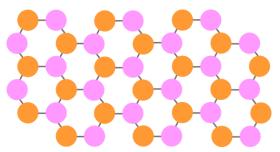


Figure 1: TMD structure MX<sub>2</sub>; M=Transition Metal (Mo) X=Chalcogen (S)

Bulk  $MoS_2$  has an indirect band gap, while in monolayer form the band gap becomes direct. The direct band gap is formed due to the removal of interlayer interactions meaning electrons are confined to a single plane. Mechanically  $MoS_2$  monolayers are highly flexible, they have a higher breaking strength than commonly used flexible plastics, and stiffness comparable to steel.<sup>1</sup>

The applications of  $MoS_2$  are numerous and wide-ranging, with interest from fields such as optoelectronics, water treatment, solar energy, and biomedicine. For the material to perform at its highest capability it must be defect free. Defects, stress, and strain can significantly influence the properties of the material. For example, removing defects responsible for non-radiative recombination allows photoluminescence quantum yields of over 95% to be realised.<sup>2</sup> Correlative Raman and photoluminescence (PL) imaging can be used to provide highly detailed sample information without any sample preparation or damage.

Using a confocal Raman microscope both Raman and PL maps can be obtained on the same sample area. In this Application Note, CVD-grown  $MoS_2$  flakes are analysed using Raman and PL imaging with the RM5 Raman Microscope.

#### Materials and Methods

An Edinburgh Instruments RM5 Raman Microscope equipped with a 532 nm laser and two gratings; 300 gr/mm for PL and 1800 gr/mm for Raman, was used to analyse MoS<sub>2</sub> flakes. The MoS<sub>2</sub> flakes were Chemical Vapor Deposition (CVD) grown on sapphire, Figure 2.

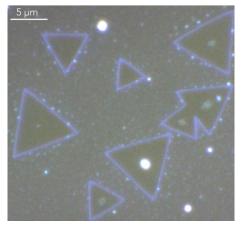


Figure 2: Reflected darkfield image of MoS<sub>2</sub> flakes.

The sample exhibits a strong Raman and PL response allowing short exposure times on the CCD camera to be used to obtain high quality spectra. During mapping When exposure times of < 0.5 seconds can be used, Fast Mapping is employed to greatly reduce mapping times. For Raman imaging an exposure time of 0.03 s was used, since PL was more intense 0.01 s exposure times could be used.

#### Raman Imaging

The Raman spectrum of  $MoS_2$  consists of two characteristic bands termed the  $E_{2g}$  band at 385 cm<sup>-1</sup> and the  $A_{1g}$  band at 405 cm<sup>-1</sup>, Figure 3. The  $E_{2g}$  band is due to in-plane vibrations, and the  $A_{1g}$  band attributed to out-of-plane vibrations.

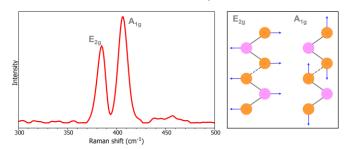


Figure 3: Raman spectrum of thin layer MoS<sub>2</sub> (left) highlighting the E<sub>2g</sub> and A<sub>1g</sub> bands and their associated vibrations (right).

# High-Resolution Raman & PL Imaging of MoS<sub>2</sub>

AN\_R15; Angela Flack



These bands are indicative of the number of layers present and the number of layers present in  $MoS_2$  crystals is critical to their function in optoelectronic devices. From Figure 4 it can be seen the  $E_{2g}$  and  $A_{1g}$  band intensity are indicative of the number of layers present, as the number of layers increases the two bands become farther apart due to inter-layer vibrations. Crystals shown in a solid single pink colour are monolayer  $MoS_2$ . While the bright spots in the centre of some of the crystals indicates the presence of multilayer  $MoS_2$ .

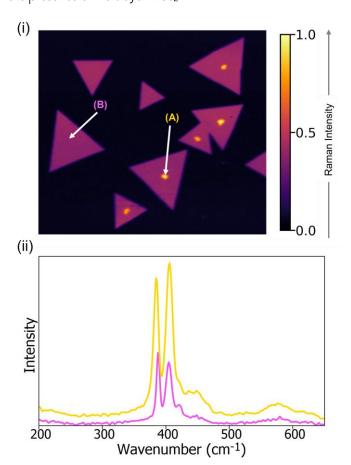


Figure 4: (i) Raman intensity imaging of MoS₂ crystals using the E₂g and A₁g bands. (ii) Raman spectra from areas of low and high Raman signal.

The splitting between the two bands was plotted, Figure 5(i). Compared to bulk  $MoS_2$  the  $E_{2g}$  band is red shifted as the crystal becomes monolayer, this shift is due to Coulomb interactions between Mo atoms which are enhanced with increased layers.<sup>3</sup> The  $A_{1g}$  band is blue shifted as the  $MoS_2$  becomes monolayer, resulting in the centre of the triangles exhibiting an increased splitting value.

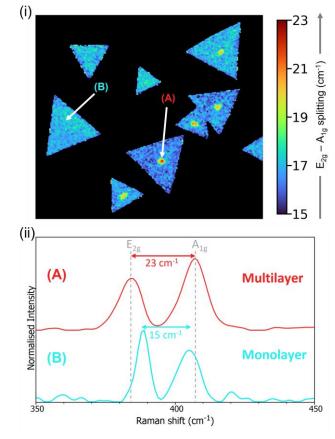


Figure 5: (i) Raman imaging displaying the band splitting of  $E_{2g}$ - $A_{1g}$ (ii) Normalised Raman spectra from monolayer and multilayer  $MoS_2$  highlighting the change in splitting.

The  $E_{2g}$  band is particularly sensitive to strain in the sample, when under strain the band is red shifted revealing the two bottom crystals highlighted in the white box are under more strain than the others, Figure 6. Raman mapping reveals these samples contain both monolayer and multilayer  $MoS_2$  whilst showing some strain.

# High-Resolution Raman & PL Imaging of MoS<sub>2</sub>

AN\_R15; Angela Flack



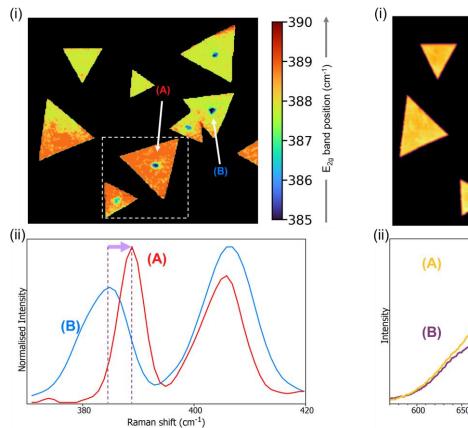


Figure 6: (i) Raman imaging displaying the band position of the  $E_2$  band and (ii) Normalised Raman spectra highlighting  $E_2$  band position.

### Photoluminescence Imaging

Excitation using a 532 nm laser offers both Raman and PL signal meaning the two phenomena can be seen in a single spectrum, therefore it is technically possible to perform the two techniques simultaneously. However, to obtain the best quality spectra different gratings are used for the two measurement types. For PL imaging a low groove density grating is used to obtain the broad PL peak. The intensity of the PL band decreases with increasing layer number and bulk  $MoS_2$  produces no PL signal at all. This effect is due to the indirect to direct band gap transition which occurs as we go from bulk  $MoS_2$  to few-layer  $MoS_2$ . The Raman and PL intensity maps should therefore be opposites of one another when comparing areas of monolayer and multilayer  $MoS_2$  which can be observed in Figure 7.

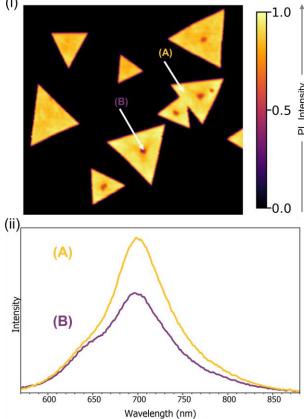


Figure 7: (i) PL intensity map of  $MoS_2$  crystals. (ii) PL spectra from an area of high and low intensity.

The PL peak is also sensitive to strain, doping, and defects, its position was plotted in Figure 8. The peak position is indicative of strain in the sample, corresponding with the  $E_{2g}$  band from Raman imaging the bottom two crystals reveal strain. Additionally, PL mapping reveals some strain in the two crystals which appear to be growing into each other, highlighted in the white box.

PL mapping offers agreement with the Raman data showing this sample has been well grown as largely monolayer crystals with little strain present.

#### **APPLICATION NOTE**

# High-Resolution Raman & PL Imaging of MoS<sub>2</sub>

AN\_R15; Angela Flack



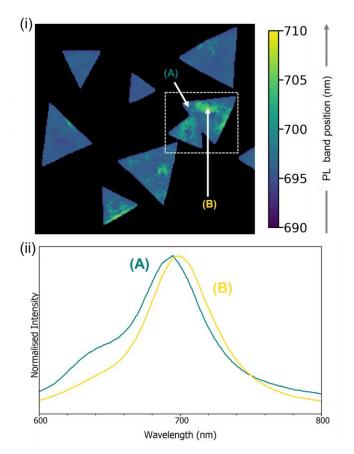


Figure 8: (i) PL imaging of MoS $_2$  crystals using peak position. (ii) Normalised PL spectra from an area of high and low intensity

## Conclusion

This application note demonstrates how the RM5 Raman Microscope can provide high resolution imaging of 2D transition metal dichalcogenides. Raman and PL imaging can be used correlatively to gain detailed information on the sample's layers, as well as indicating strain, defects, and doping.

### References

- 1. Pei, J. et al. Exciton and Trion Dynamics in Bilayer MoS2. Small 11, 6384–6390 (2015).
- 2. Amani, M. et al. Near-unity photoluminescence quantum yield in MoS2. *Science (80-. ).* **350**, 1065–1068 (2015).
- Golovynskyi, S. et al. Exciton and trion in few-layer MoS2: Thickness- and temperature-dependent photoluminescence. Appl. Surf. Sci. 515, (2020).



For more information, please contact:

+44 (0) 1506 425 300 sales@edinst.com www.edinst.com