APPLICATION NOTE

Transient Absorption and Singlet Oxygen Phosphorescence of Rose Bengal Photosensitiser

AN_TA01; Maria Tesa



Introduction

A photosensitiser (PS) is a chemical species that initiates a reaction upon absorption of light. Photosensitisers vary widely in their chemical composition and practical applications, but they all follow the same mechanism: the PS absorbs light and reaches a triplet excited state. This triplet state, which has a typical lifetime of microseconds, interacts with a target species, transferring energy that induces further reactions. The PS returns to its ground state and can take part in a new photochemical activation.

A widespread application of PSs is the generation of singlet oxygen ($^{1}O_{2}$) in photodynamic therapy, a treatment for some types of cancer and skin conditions. The PS is introduced to the patient's body and excited with light into a singlet state, which transitions into a triplet state via intersystem crossing (Figure 1). Energy transfer then takes place between the excited PS and ground-state molecular oxygen ($^{3}O_{2}$) to generate $^{1}O_{2}$. Singlet oxygen is extremely reactive and destroys the target cells, either by direct reaction or by generating oxidised products that react further. The process can be probed by different types of spectroscopy: fluorescence or phosphorescence of the photosensitiser, transient absorption (TA) of the triplet state, and $^{1}O_{2}$ phosphorescence.

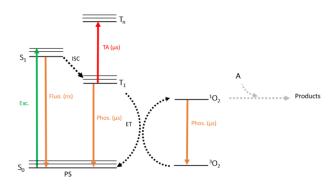


Figure 1 Jablonski diagram showing excitation of a photosensitiser and singlet oxygen generation. The PS first absorbs light (Exc.) into S_1 , then undergoes intersystem crossing (ISC) into T_1 . T_1 transfers energy (ET) to 3O_2 to generate 1O_2 , which reacts with other species A to generate photodynamic therapy products.

There is extensive research to find suitable PSs for photodynamic therapy that produce high concentrations of ${}^{1}O_{2}$. The ideal PS molecule yields a high T₁ state population upon excitation. In addition, the T₁ state should be long-lived and not quenched by species other than ${}^{3}O_{2}$. Optimising the PS requires understanding its triplet formation mechanism, which usually involves characterising the spectrum and lifetime of the T₁ state. This is achieved via nanosecond transient absorption (TA), which probes the absorption of T₁ as it is excited into higher-lying triplet states (Figure 1). In addition, direct detection of ${}^{1}O_{2}$ is often employed to characterise the yield of energy transfer from T₁ to ${}^{3}O_{2}$.

In this application note, an Edinburgh Instruments LP980 Transient Absorption Spectrometer is employed to study the triplet state of rose bengal, a common ${}^{1}O_{2}$ photosensitiser. The same instrument is also used to directly detect ${}^{1}O_{2}$ by laser-induced phosphorescence in the NIR range, eliminating the need for a separate photoluminescence spectrometer.

Materials and methods

A solution was prepared by dissolving the photosensitiser rose bengal in acetonitrile to an optical density of 0.1 at 510 nm and ~0.5 at 560 nm and loaded into a quartz degassing cuvette. The solution was either saturated with air or degassed by three consecutive freezepump-thaw cycles for oxygen-free measurements.

An Edinburgh Instruments LP980 Transient Absorption Spectrometer (Figure 2) was used to study the time-resolved absorption and luminescence of the PS in the visible range. A frequency-doubled Nd:YAG laser (Litron Nano S 130-10) provided 532 nm excitation pulses of 5-7 ns with a frequency of 10 Hz. The standard pulsed Xe lamp in the LP980 was used as the probe. Kinetic and spectral data were acquired with a PMT-900 and intensified CCD (ICCD) detector, respectively.

Time-resolved ${}^{1}O_{2}$ phosphorescence was studied in the LP980 with the same pump laser, replacing the visible PMT-900 with an NIR-sensitive PMT-1400 and photon-counting Multichannel Scaling (MCS) electronics.



Figure 2 LP980 Transient Absorption Spectrometer.

TA and Fluorescence Spectra of Rose Bengal

Figure 3 presents TA and fluorescence spectra from rose bengal in acetonitrile. TA spectra measure the absorption from the transient triplet state, while laser-induced fluorescence (LIF) spectra measure photoluminescence emission. Data were acquired for a degassed or air-saturated solution. The presence of oxygen in the air-saturated sample shortens the lifetime of the triplet, which is quenched into its ground state, generating ${}^{1}O_{2}$ at the same time. Such strong

APPLICATION NOTE

Transient Absorption and Singlet Oxygen Phosphorescence of Rose Bengal Photosensitiser



AN_TA01; Maria Tesa

dependence of the triplet lifetime on air concentration is characteristic of an efficient PS.

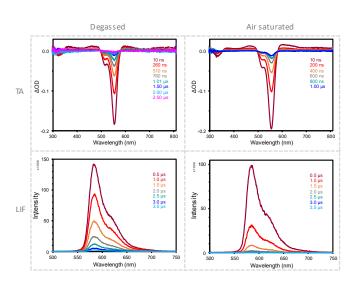


Figure 3 Transient absorption (TA) and fluorescence (LIF) spectra from rose bengal in acetonitrile acquired in an LP980 spectrometer under degassed (left) and air-saturated (right) conditions. Measurement parameters: $\lambda_{pump} = 532 \text{ nm}$, $E_{pump} = 10 \text{ mJ/pulse}$, $\Delta \lambda_{probe} = 1 \text{ nm}$, repetition rate = 10 Hz. The delays between ICCD acquisition and pump laser are indicated in the plots.

Without prior knowledge of the PS, the LIF spectra in Figure 3 could be interpreted as phosphorescence from the T_1 state. However, the results do not match the phosphorescence spectrum of rose bengal,¹ and are identical to its fluorescence spectrum. The luminescence is too long-lived to be prompt fluorescence from the S_1 state; besides this, it is quenched by oxygen, suggesting that it arises from a triplet. A potential explanation is delayed fluorescence arising from either triplet-triplet annihilation or reverse intersystem crossing² upon interaction with ¹O₂.

TA Decays of Rose Bengal

The triplet relaxation dynamics can be studied in detail with timeresolved TA traces. Figure 4 presents examples using the PMT detector in LP980. Both the negative ground state bleach (a) and the positive TA (b) components of the spectrum are longer lived in the absence of oxygen, confirming the qualitative trend observed in Figure 3.

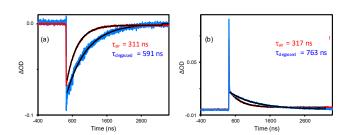


Figure 4 Transient absorption decays from rose bengal in acetonitrile acquired in LP980 spectrometer at (a) 520 nm and (b) 440 nm under degassed (blue) and air-saturated (red) conditions. Measurement parameters: $\lambda_{pump} = 532 \text{ nm}$, $E_{pump} = 10 \text{ mJ/pulse}$, $\Delta\lambda_{probe} = 0.70 \text{ nm}$, repetition rate = 10 Hz. Fit results to a single exponential decay model are shown in the figure.

NIR Phosphorescence Decays of ¹O₂

Transient absorption of the triplet state provides an insight into the photophysics of the process, and it can be used to screen multiple PS candidates. However, direct detection of singlet oxygen phosphorescence in the NIR remains the best tool for evaluating the PS performance. Such experiments are typically performed in a dedicated photoluminescence spectrometer like the Edinburgh Instruments FLS1000.³ The LP980 removes the need for a separate spectrometer as it can be fitted with a photon-counting NIR PMT and electronics identical to those in FLS1000. This is not only convenient for the user, but it also allows comparison between the TA and NIR phosphorescence datasets with less uncertainty: the same pump laser, sample chamber, and detection monochromator are used so the two experiments are more directly comparable.

A ${}^{1}O_{2}$ phosphorescence decay acquired in the LP980 is shown in Figure 5. The data was acquired in a few minutes thanks to the high sensitivity of the photon-counting NIR PMT detector. The decay fits to a single exponential model with a lifetime of 67 microseconds.

APPLICATION NOTE

Transient Absorption and Singlet Oxygen Phosphorescence of Rose Bengal Photosensitiser



AN_TA01; Maria Tesa

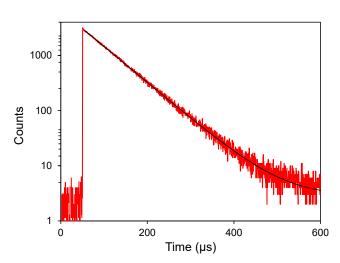


Figure 5 NIR phosphorescence decay from ${}^{1}O_{2}$ acquired in LP980 exciting with 532 nm Nd:YAG laser pulses (laser scatter subtracted from the data) and multichannel scaling detection. The exponential fit result is shown in the graph. Measurement conditions: laser energy = 5 nJ/pulse, repetition rate = 10 Hz, λ_{em} = 1275 nm, $\Delta\lambda_{em}$ = 3 nm, 20 minutes acquisition.

Conclusion

The LP980 Transient Absorption Spectrometer was successfully used to acquire the spectrum and lifetime of the rose bengal triplet state, as well as visible photoluminescence from rose bengal and NIR phosphorescence from its singlet oxygen product. A complete photophysical characterisation of the photosensitiser and singlet oxygen is possible thanks to the versatile configuration options of the LP980.

References

1. A. Penzkofer et al., Room temperature phosphorescence lifetime and quantum yield of erythrosine B and rose bengal in aerobic alkaline aqueous solution, J. Lumin. 132 1055-1062 (2012)

For more information, please contact:



For more information, please contact:

+44 (0) 1506 425 300 sales@edinst.com www.edinst.com

- M. Scholz et al., Singlet oxygen-sensitized delayed fluorescence of common water-soluble photosensitizers, Photochem. Photobiol. Sci. 12 1873-1884 (2013)
- Tesa, Maria. Detection of Singlet Oxygen by Photoluminescence Spectroscopy, Edinburgh Instrument Application Note AN_P85 (2023)